

Xeno-Free, Viral Inactivated, Exosome-Depleted for EV Production

AventaCell BioMedical Corp. has developed an exosome depletion process to remove human platelet lysate (hPL)-derived exosomes. Exosome-Depleted UltraGRO™-PURE GI (ED UG-P GI) is able to support human MSC cell viability to secrete abundant extracellular vesicles (EVs) without compromising phenotype over the culture period. Moreover, gamma irradiation processing of the product is used as a pathogen reduction treatment (PRT) for viral inactivation, to comply with regulatory guidance for clinical research and development.

Benefits of Exosome-Depleted UltraGRO™-PURE GI

- Xeno-free with >95% nanoparticle removal from the hPL supplement
- Minimal hPL nanoparticle contamination
- MSCs cultured with the depleted supplement remain highly viable with stable phenotype markers throughout the culture period
- GMP Exosome-Depleted UltraGRO™-PURE GI to produce clinical grade hMSC-derived EVs
- Gamma irradiation processing is accepted by regulatory agencies as a validated PRT

High depletion rate was performed in the manufacturing process

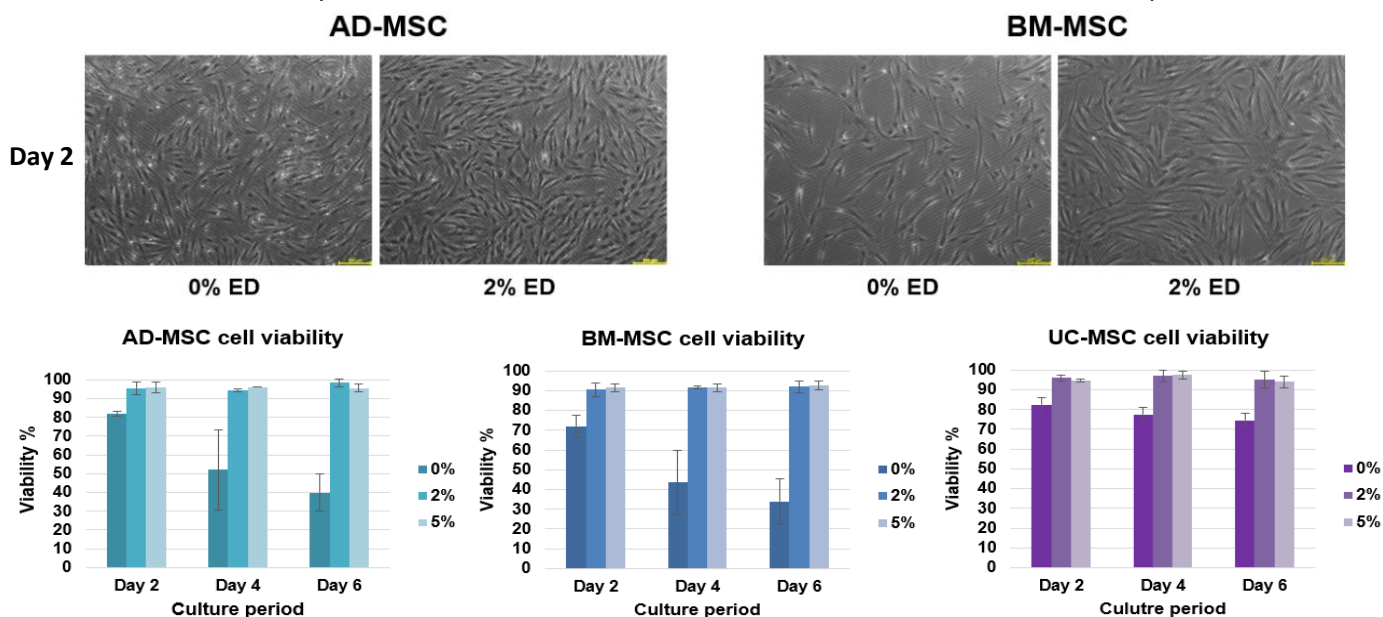
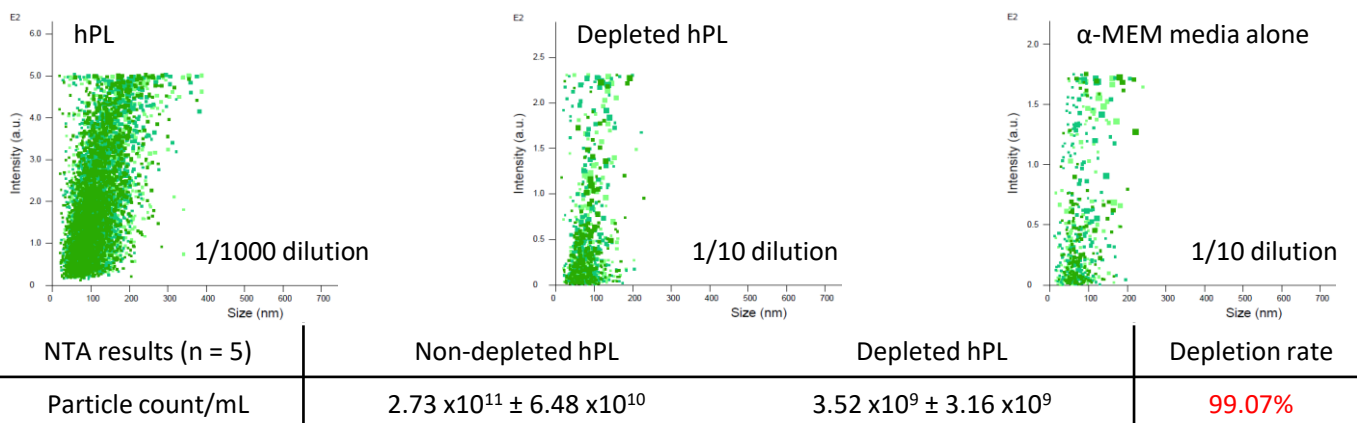


Fig 1. Highly depletion rate was performed, and ED UG-P GI is able to support MSCs for long-term EV production.

Particle in medium	α -MEM media alone	2% ED UG-P GI	5% ED UG-P GI	<i>Minimal interference</i>
Particle count/mL	$1.93 \times 10^8 \pm 1.27 \times 10^7$	$3.51 \times 10^8 \pm 1.31 \times 10^7$	$4.32 \times 10^8 \pm 5.54 \times 10^7$	

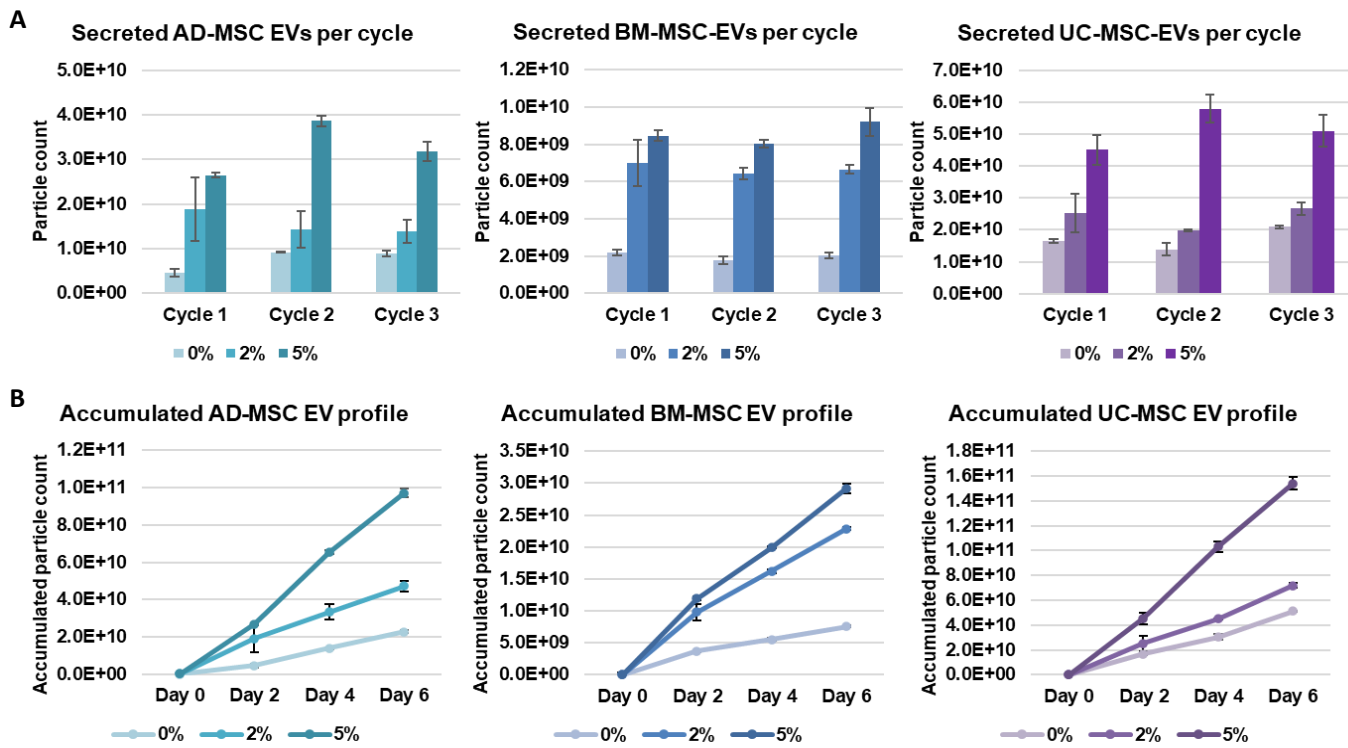


Fig.2: After MSCs reached 50-70% confluency in a T75 flask, the culture media was changed to ED UG-P GI supplemented culture media, and the media refresh was performed every 2 days as one culture cycle. (A) MSC-derived EVs secretion and their (B) accumulation profile were measured by NTA. (C) The MSC specific phenotype was not altered throughout the culture period up to 14 days.

C MSC phenotype characterization

Culture period	CD73	CD90	CD105	CD34
Day 10	97.4%	99.4%	99.7%	1.2%
Day 14	97.9%	99.8%	97.6%	1.8%

Billions of MSC-derived EVs can be easily acquired for further applications!



Specifications	Acceptance
Appearance	Sligh yellow
Mycoplasma	Negative
Endotoxin	< 10 EU/mL
Sterility	No growth
pH	6.5 – 8.5
Osmolarity	270 – 330 mOsm/kg
Particle depletion rate	> 95%
Cell assay	Support MSC culture
Gamma irradiation dose	25 – 40 kGy

Ordering Information

Product Number	Product	Bottle Size (mL)
HPCHEFRLI05	Exosome-Depleted UltraGRO™-PURE GI	50
HPCHEFRLI10		100
HPCHEFRLI50		500
HPCHEFGLI05	Exosome-Depleted UltraGRO™-PURE GI (GMP grade)	50
HPCHEFGLI10		100
HPCHEFGLI50		500

CliniSciences Group

Austria

Company: CliniSciences GmbH
Address: Sternwartestrasse 76, A-1180
Wien - Austria
Telephone: +43 720 115 580
Fax: +43 720 115 577
Email: oesterreich@clinisciences.com
Web: <https://www.clinisciences.com>



Belgium

Company: CliniSciences S.R.L
Address: Avenue Stalingrad 52, 1000
Brussels - Belgium
Telephone: +32 2 31 50 800
Fax: +32 2 31 50 801
Email: belgium@clinisciences.com
Web: <https://www.clinisciences.com>



Denmark

Company: CliniSciences ApS
Address: Oesterbrogade 226, st. 1,
Copenhagen, 2100 - Denmark
Telephone: +45 89 888 349
Fax: +45 89 884 064
Email: danmark@clinisciences.com
Web: <https://www.clinisciences.com>



Finland

Company: CliniSciences ApS
Address: Oesterbrogade 226, st. 1,
Copenhagen, 2100 - Denmark
Telephone: +45 89 888 349
Fax: +45 89 884 064
Email: suomi@clinisciences.com
Web: <https://www.clinisciences.com>



France

Company: CliniSciences S.A.S
Address: 74 Rue des Suisses, 92000
Nanterre - France
Telephone: +33 9 77 40 09 09
Fax: +33 9 77 40 10 11
Email: info@clinisciences.com
Web: <https://www.clinisciences.com>



Germany

Company: Biotrend Chemikalien GmbH
Address: Wilhelm-Mauser-Str. 41-43,
50827 Köln - Germany
Telephone: +49 221 9498 320
Fax: +49 221 9498 325
Email: info@biotrend.com
Web: <https://www.biotrend.com>



Iceland

Company: CliniSciences ApS
Address: Oesterbrogade 226, st. 1,
Copenhagen, 2100 - Denmark
Telephone: +45 89 888 349
Fax: +45 89 884 064
Email: island@clinisciences.com
Web: <https://www.clinisciences.com>



Ireland

Company: CliniSciences Limited
Address: Ground Floor, 71 lower Baggot street
Dublin D02 P593 - Ireland
Telephone: +353 1 6971 146
Fax: +353 1 6971 147
Email: ireland@clinisciences.com
Web: <https://www.clinisciences.com>



Italy

Company: CliniSciences S.r.l
Address: Via Maremmana inferiore 378
Roma 00012 Guidonia Montecelio - Italy
Telephone: +39 06 94 80 56 71
Fax: +39 06 94 80 00 21
Email: italia@clinisciences.com
Web: <https://www.clinisciences.com>



Netherlands

Company: CliniSciences B.V.
Address: Kraaijenhoffstraat 137A,
1018RG Amsterdam, Netherlands
Telephone: +31 85 2082 351
Fax: +31 85 2082 353
Email: nederland@clinisciences.com
Web: <https://www.clinisciences.com>



Norway

Company: CliniSciences ApS
Address: Oesterbrogade 226, st. 1,
Copenhagen, 2100 - Denmark
Telephone: +45 89 888 349
Fax: +45 89 884 064
Email: norge@clinisciences.com
Web: <https://www.clinisciences.com>



Poland

Company: CliniSciences sp.Z.o.o.
Address: ul. Rotmistrza Witolda Pileckiego 67
lok. 200 - 02-781 Warszawa -Poland
Telephone: +48 22 307 0535
Fax: +48 22 307 0532
Email: polska@clinisciences.com
Web: <https://www.clinisciences.com>



Portugal

Company: Quimigen Unipessoal LDA
Address: Rua Almada Negreiros, Lote 5, Loja 14,
2615-275 Alverca Do Ribatejo - Portugal
Telephone: +351 30 8808 050
Fax: +351 30 8808 052
Email: info@quimigen.com
Web: <https://www.quimigen.pt>



Spain

Company: CliniSciences Lab Solutions
Address: C/ Hermanos del Moral 13
(Bajo E), 28019, Madrid - Spain
Telephone: +34 91 269 40 65
Fax: +34 91 269 40 74
Email: espana@clinisciences.com
Web: <https://www.clinisciences.com>



Sweden

Company: CliniSciences ApS
Address: Oesterbrogade 226, st. 1,
Copenhagen, 2100 - Denmark
Telephone: +45 89 888 349
Fax: +45 89 884 064
Email: sverige@clinisciences.com
Web: <https://www.clinisciences.com>



Switzerland

Company: CliniSciences Limited
Address: Marktgasse 18 8302 Kloten -
Switzerland
Telephone: +41 (044) 805 76 81
Fax: +41 (044) 805 76 75
Email: switzerland@clinisciences.com
Web: <https://www.clinisciences.com>



UK

Company: CliniSciences Limited
Address: 11 Progress Business center, Whittle
Parkway, SL1 6DQ Slough- United Kingdom
Telephone: +44 (0)1753 866 511
or +44 (0) 330 684 0982
Fax: +44 (0)1753 208 899
Email: uk@clinisciences.com
Web: <https://www.clinisciences.com>



USA

Company: Biotrend Chemicals LLC
Address: c/o Carr Riggs Ingram,
500 Grand Boulevard, Suite 210 Miramar
Beach, FL 32550- USA
Telephone: +1 850 650 7790
Fax: +1 850 650 4383
Email: info@biotrend-usa.com
Web: <https://www.biotrend-usa.com>

