



READYTECTOR.
easy, quick and clear

Distributed by:

CliniSciences Group

NEW

The all-in-one Solution for Western blotting!

The all-in-one Solution for Western blotting!

ReadyTector® is the all-in-one detection solution for Western blotting.

ReadyTector® contains everything you need for fast, all-in-one immunodetection. Only a specific primary antibody has to be added. All-in-one means that one solution contains everything you need, and the entire process takes place in a single work step – blocking and binding of the primary and secondary antibody occur simultaneously. The next step is just a washing sequence using the special ReadyTector® Wash Buffer.

The ReadyTector® solution with your valuable primary antibody can be re-used up to 5 times.


ReadyTector® kits include the all-in-one solution and 10× Wash Buffer, each in 40 mL, 120 mL and 500 mL bottles.

For detection with chemiluminescence we recommend the specific ReadyTector® Chemiluminescent Substrate.

Despite the fast, all-in-one incubation process, ReadyTector® reduces background, allowing users to generate clear, distinct bands suitable for publishing.

**ReadyTector® –
easy, quick and clear**

Comparable results obtained with standard Western blotting protocol or ReadyTector® protocol and ReadyTector® Chemiluminescent Substrate

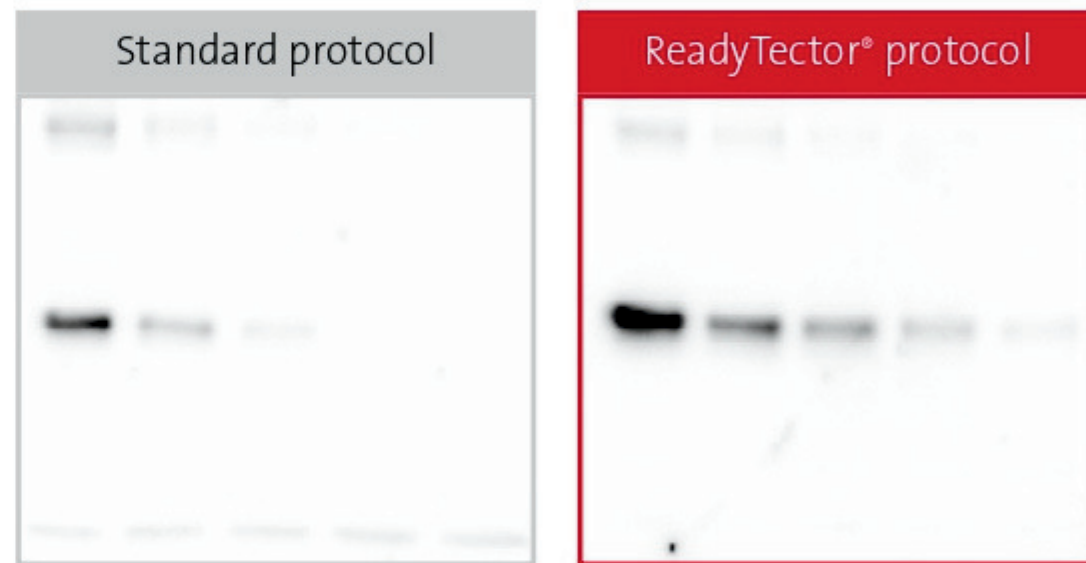

Lanes 1–5 contain 22, 11, 5.5, 2.8, 1.4 ng Alpha-1-Antitrypsin spiked in cell lysates, respectively.

Protein detection: Mouse anti-A₁AT Clone 1AT (Biotrend), 0.2 µg / mL on nitrocellulose.

The primary antibody for A₁AT should only show one band. Additional weak signals show non-specific extra bands.

Distributed by:

CliniSciences Group

Easy, quick and clear

Easy – It is very easy to switch from the classic immunodetection procedure with several incubation steps to the faster and more convenient ReadyTector® method.

With ReadyTector®, immunodetection will be accelerated and simplified. All other working

steps and materials will remain unchanged. You can also use the same primary antibodies. Blotting procedure, membranes and signal detection do not have to be changed.

No new device or equipment is necessary. ReadyTector® is easy to use and universal.

Quick – Fewer working and washing steps save you time and provide results more quickly.

The ReadyTector® solution contains the blocker and secondary antibodies and thus it can be used directly without any dilution or pre-incubation steps. You simply add your specific primary antibody.

Clear – ReadyTector® reduces background and enables clear and sharp bands for most applications. The results are suitable for publishing.

ReadyTector® protocol



ReadyTector® saves several washing and incubation steps. Therefore a time saving of several hours can be obtained.

Run your gel and blot and detect the membrane in a single morning. You get the same great results with ReadyTector® – simply much faster.

Standard protocol



Distributed by:

CliniSciences Group



CliniSciences Group

Austria

Company: CliniSciences GmbH
 Address: Sternwartestrasse 76, A-1180
 Wien - Austria
 Telephone: +43 720 115 580
 Fax: +43 720 115 577
 Email: osterreich@clinisciences.com
 Web: <https://www.clinisciences.com>



Belgium

Company: CliniSciences S.R.L.
 Address: Avenue Stalingrad 52, 1000
 Brussels - Belgium
 Telephone: +32 2 31 50 800
 Fax: +32 2 31 50 801
 Email: belgium@clinisciences.com
 Web: <https://www.clinisciences.com>



Denmark

Company: CliniSciences ApS
 Address: Oesterbrogade 226, st. 1,
 Copenhagen, 2100 - Denmark
 Telephone: +45 89 888 349
 Fax: +45 89 884 064
 Email: danmark@clinisciences.com
 Web: <https://www.clinisciences.com>



Finland

Company: CliniSciences ApS
 Address: Oesterbrogade 226, st. 1,
 Copenhagen, 2100 - Denmark
 Telephone: +45 89 888 349
 Fax: +45 89 884 064
 Email: suomi@clinisciences.com
 Web: <https://www.clinisciences.com>



France

Company: CliniSciences S.A.S
 Address: 74 Rue des Suisses, 92000
 Nanterre - France
 Telephone: +33 9 77 40 09 09
 Fax: +33 9 77 40 10 11
 Email: info@clinisciences.com
 Web: <https://www.clinisciences.com>



Germany

Company: Biotrend Chemikalien GmbH
 Address: Wilhelm-Mauser-Str. 41-43,
 50827 Köln - Germany
 Telephone: +49 221 9498 320
 Fax: +49 221 9498 325
 Email: info@biotrend.com
 Web: <https://www.biotrend.com>



Iceland

Company: CliniSciences ApS
 Address: Oesterbrogade 226, st. 1,
 Copenhagen, 2100 - Denmark
 Telephone: +45 89 888 349
 Fax: +45 89 884 064
 Email: island@clinisciences.com
 Web: <https://www.clinisciences.com>



Ireland

Company: CliniSciences Limited
 Address: Ground Floor, 71 lower Baggot street
 Dublin D02 P593 - Ireland
 Telephone: +353 1 6971 146
 Fax: +353 1 6971 147
 Email: ireland@clinisciences.com
 Web: <https://www.clinisciences.com>



Italy

Company: CliniSciences S.r.l
 Address: Via Maremmana inferiore 378
 Roma 00012 Guidonia Montecelio - Italy
 Telephone: +39 06 94 80 56 71
 Fax: +39 06 94 80 00 21
 Email: italia@clinisciences.com
 Web: <https://www.clinisciences.com>



Netherlands

Company: CliniSciences B.V.
 Address: Gaetano Martinolaan 85, 6229
 GS Maastricht - Netherlands
 Telephone: +31 85 2082 351
 Fax: +31 85 2082 353
 Email: nederland@clinisciences.com
 Web: <https://www.clinisciences.com>



Norway

Company: CliniSciences ApS
 Address: Oesterbrogade 226, st. 1,
 Copenhagen, 2100 - Denmark
 Telephone: +45 89 888 349
 Fax: +45 89 884 064
 Email: norge@clinisciences.com
 Web: <https://www.clinisciences.com>



Poland

Company: CliniSciences sp. z o.o.
 Address: ul. Rotmistrza Witolda Pileckiego 67
 lok. 200 - 02-781 Warszawa - Poland
 Telephone: +48 22 307 0535
 Fax: +48 22 307 0532
 Email: polaka@clinisciences.com
 Web: <https://www.clinisciences.com>



Portugal

Company: Quimigen Unipessoal LDA
 Address: Rua Almada Negreiros, Lote 5, Loja 14,
 2615-275 Alverca Do Ribatejo - Portugal
 Telephone: +351 30 8808 050
 Fax: +351 30 8808 052
 Email: info@quimigen.com
 Web: <https://www.quimigen.pt>



Spain

Company: CliniSciences Lab Solutions
 Address: C/ Hermanos del Moral 13
 (Bajo E), 28019, Madrid - Spain
 Telephone: +34 91 269 40 65
 Fax: +34 91 269 40 74
 Email: espana@clinisciences.com
 Web: <https://www.clinisciences.com>



Sweden

Company: CliniSciences ApS
 Address: Oesterbrogade 226, st. 1,
 Copenhagen, 2100 - Denmark
 Telephone: +45 89 888 349
 Fax: +45 89 884 064
 Email: sverige@clinisciences.com
 Web: <https://www.clinisciences.com>



Switzerland

Company: CliniSciences Limited
 Address: Marktgasse 18 8302 Kloten -
 Switzerland
 Telephone: +41 (044) 805 76 81
 Fax: +41 (044) 805 76 75
 Email: switzerland@clinisciences.com
 Web: <https://www.clinisciences.com>



UK

Company: CliniSciences Limited
 Address: 11 Progress Business center, Whittle
 Parkway, SL1 6DQ Slough - United Kingdom
 Telephone: +44 (0)1753 866 511
 or +44 (0) 330 684 0982
 Fax: +44 (0)1753 208 899
 Email: uk@clinisciences.com
 Web: <https://www.clinisciences.com>



USA

Company: Biotrend Chemicals LLC
 Address: c/o Carr Riggs Ingram,
 500 Grand Boulevard, Suite 210 Miramar
 Beach, FL 32550- USA
 Telephone: +1 850 650 7790
 Fax: +1 850 650 4383
 Email: info@biotrend-usa.com
 Web: <https://www.biotrend-usa.com>

